Manutan

By your side, every day

A227360

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COMBINED SHOWER AND EYEBATH

TECHNICAL SPECIFICATION

	A227360
Type of control	Main
Shower head material	Steel
Base material	Stainless steel
Frame material	Stainless steel
Head Ø	255 mm
Base Ø	145 mm
O/A height	2450 mm

- Head: 304 stainless steel shower head.
- Eye Wash Nozzle: green ABS spray, others are 304 stainless steel with 10" waste water recycle bowl
- Shower Valve: 1" 304 stainless steel ball valve
- Eye Wash Valve: 1/2" 304 stainless steel ball valve
- Supply: 1 1/4" FNPT
- Waste: 1 1/4" FNPT

PRODUCT CONCEPT

Emergency eye wash and shower units are designed to rinse contaminants from the user's eyes, face or body. As such, these units are forms of first aid equipment to be used in the event of an accident. However, they are not a substitute for primary protective devices (including eye and face protection and protective clothing) or for safety procedures when handling hazardous materials.

A227360 Stainless Steel Combination Eye Wash/Shower main material is 304 stainless steel, which is particularly suitable for use with weak acids and alkalis. If strong acids, alkalis or chlorides are present in the environment then we recommend using the 316 stainless steel or ABS eye wash.

LOCATION

- In general, the ANSI standard requires that emergency equipment be installed within 10 seconds walking distance from the location of a hazard (approximately 55 feet).
- The equipment must be installed on the same level as the hazard (i.e. accessing the equipment should not require going up or down stairs or ramps).
- The path of travel from the hazard to the equipment should be free of obstructions and as direct as possible.
- The location of the emergency equipment must be marked with a highly visible sign.

ORIGINAL WATER

Drinking water or filtered water

HYDRAULIC PRESSURE AND FLOW RATE

If the pressure exceeds 0.2-0.6MPa, please install the relief valve to protect the user's eyes. Eye wash flow $\geq 11.4\text{L/Min}$, shower flow $\geq 75.7\text{L/Min}$.

TRAINING

Simply installing emergency equipment is not sufficient means of ensuring worker safety. It is also very important that employees are trained in the location and proper use of emergency equipment. Research shows that after an incident has occurred, rinsing eyes within the first ten seconds is essential. Therefore, employees at the highest risk of damaging their eyes in each department must be trained regularly. All employees must know the location of the emergency equipment and be aware that quick and effective rinsing is important in an emergency.

Eye/Face Wash

- The sooner the injured employee's eyes are rinsed, the lower the risk of damage. Every second is important when preventing permanent damage to save time for medical treatment.
- All employees must be reminded that this equipment is only to be used in emergencies, tampering with the equipment may cause a malfunction.
- In emergencies, the afflicted may be unable to open their eyes. Employees may feel pain, anxiousness and loss. They may need others' help to reach the equipment and use it.
- Push the handle to spray the liquid.
- When liquid sprays, put the injured employee's left hand on the left nozzle, and the right hand on the right nozzle.
- Put the injured employee's head over the eye wash bowl which is hand controlled,
- When rinsing the eyes, use both hands' thumb and index finger to open the eyelids, rinsing for at least 15 minutes.
- After rinsing, seek medical treatment immediately
- Security and supervisory personnel must be notified that the equipment has been used.

Shower

- Use the pull rod to start the liquid flow.
- The injured should stand in the water flow once it has started.
- Ensure affected areas are in the water flow.
- Do not rinse by hand, to avoid further injury.

Note: If chemicals which react dangerously with water are present, an alternative harmless liquid will be provided. Special eye drops should also be used.

INSTRUCTIONS

► Figure 1 of this product can be found at the end of this document.

1	Optional Supply: 1 1/4" FNPT
2	10" 304 stainless steel shower head
3	Shower Valve: 1" 304 stainless steel ball valve
4	Supply: 1 1/4" FNPT
5	Main Pipe: 304 stainless steel
6	Eye Wash Nozzle: green ABS spray, others are 304 stainless steel with 10" waste water recycle bowl
7	3/8" outside wire
8	Eye Wash Valve: 1/2" 304 stainless steel ball valve
9	Waste: 1 1/4" FNPT

INSTALLATION

► Figure 2 of this product can be found at the end of this document.

1	Take out the base and fix to the floor.
2	Connect the T-type drainage pipe to the base.
3	Connect the main pipe to the T-type drainage pipe.
4	Connect the main body to the main pipe, facing the user.
5	Screw the eye wash bowl onto the main body.
6	Screw the eye wash nozzle onto the main body.
7	Install the lever onto the main body valve and fasten it securely.
8	Connect the main pipe to the main body (warning tape).
9	Connect the inlet T-type to the main pipe.
10	Connect the main pipe to the inlet T-type.
11	Connect the T-type ball valve to the main pipe.
12	Connect the horizontal pipe to the T-type ball valve.
13	Connect the elbow pipe to the horizontal pipe, pointing downwards.
14	Connect the shower head to the elbow pipe.
15	Connect the pull rod and valve handle to the T-type ball valve.

Attention: Wrap every joint between pipes with PTFE tape.

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